

LLF2 Combination Cam Lock Instructions



LLF2 combination lock installed

I. Manufacturer's combination

The manufacturer has set the combination code at 0-0-0-0.

II. Resetting the combination

- Make sure the digit wheels are set at the correct code so that the housing turns freely. The cam should be engaged so that the doors of the appliance cannot open.
- Use a screwdriver or other tool to push down and hold the reset button until you have entered your new code. (Fig. 1)
- Turn the four digit wheels to your new combination code, then release the reset button.
- Make sure the housing is able to turn freely at the new setting. Code setting is complete.

III. To lock

While the lock is in the unlocked position (i.e., the housing can be turned by hand), randomly set the digit wheels to lock. (Fig. 2)

IV. To unlock

Turn the digit wheels to the correct code to unlock. (Fig. 3)

V. Use key to unlock

You can use the key to unlock if you forget the combination.

- Insert key into the barrel. (Fig. 4)
- Turn the key 90 degrees to unlock. If the key is used to unlock, the cam will also turn 90 degrees when the key is rotated. However, the digit wheels will not rotate.

