INSTRUCTIONS

Power Failure Alarm Cord, Model PF1

This ACCUCOLD power-failure line cord has a built-in power failure sensor with an 80dB loud audible alarm. It is intended for monitoring essential 120 volt AC powered cord-connected appliances for power failure events, accidental cord disconnection and circuit breaker failure. The loud signal sounds for up to 90 minutes, advising the user to reconnect power to the device.

Key Components of the Power Failure Alarm Cord

- ✓ 15 Amp hospital grade grounded three-wire plug
- ✓ Alarm housing with rechargeable battery
- ✓ AC Power ON LED indicator
- ✓ Grounded three-wire 15 Amp connector
- ✓ Built-in switch that will silence the alarm when operated in the “alarm off” position

INSTALLATION

- i. Attach the female connector to the plug of the appliance to be monitored.
- ii. Plug the cord into any live 120 volt AC outlet.

Note: If the appliance plug is disconnected from the cord, the alarm will not sound. It will sound only if power fails to reach the cord.

SAVE THIS MANUAL FOR FUTURE REFERENCE

TESTING AND OPERATION

After the cord has been plugged into an AC outlet for 24 hours to obtain a full charge, simply disconnect the cord from the AC outlet. The alarm should sound. Make certain that the switch is in the “alarm on” position.

SPECIFICATIONS

- Voltage and current: 120 VAC maximum, 15 amps maximum
- Frequency: 60 Hz
- For indoor use only
- Operating temperature range: 5°C to 40°C (41°F to 104°F)

Limited Warranty

SUMMIT products have a limited warranty period of 1 year against defects in materials and workmanship from the date of purchase. Accessory items and sensors have a limited warranty of 3 months. Repair services have a limited warranty period of 3 months against defects in materials and workmanship. SUMMIT shall, at its option, either repair or replace hardware products that prove to be defective, if a notice to that effect is received within the warranty period. SUMMIT makes no other warranties or representations of any kind whatsoever, expressed or implied, except that of title, and all implied warranties including any warranty of merchantability and fitness for a particular purpose are hereby disclaimed.

Summit, Div of Felix Storch, Inc. • Bronx, NY 10474 • Tel: 718-893-3900 • Fax: 844-478-8799

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Summit, Div of Felix Storch, Inc. • Bronx, NY 10474 • Téléphone: 718-893-3900 • Fax: 844-478-8799